

DEPARTMENT OF HEALTH CARE ACCESS AND INFORMATION FACILITIES DEVELOPMENT DIVISION

OFFICE USE ONLY APPLICATION FOR HCAI PREAPPROVAL OF **MANUFACTURER'S CERTIFICATION (OPM)** APPLICATION #: OPM-0219 **HCAI Preapproval of Manufacturer's Certification (OPM)** Renewal/Update New Type: **Manufacturer Information** Manufacturer: Belimed Infection Control Manufacturer's Technical Representative: John Nies Mailing Address: 8351 Palmetto Commerce Parkway Suite 101, Ladson, SC 29456 Telephone: (800) 451-4118 Email: john.nies@belimed.com **Product Information** Product Name: Ultrasonic Washers Product Type: Laboratory Equipment Product Model Number: 200c & UW565 General Description: Floor mounted ultrasonic washers with seismic mounting brackets **Applicant Information** Applicant Company Name: EASE LLC. Contact Person: Tiffany Tonn

"A healthier California where all receive equitable, affordable, and quality health care"

Mailing Address: 1515 FAIRVIEW AVE, STE 205, MISSOULA, MT 59801

Man

Telephone: (406) 541-3273

Title: Office Manager

Email: tiffany@easeco.com



DEPARTMENT OF HEALTH CARE ACCESS AND INFORMATION FACILITIES DEVELOPMENT DIVISION

Registered Design Professonal Preparing Engineering Recommendations
Company Name: EASE LLC
Name: Jonathan Roberson California License Number: S4197
Mailing Address: 5877 Pine Ave., Suite 210, Chino Hills, CA 91709
Telephone: (951) 295-1892
HCAI Special Seismic Certification Preapproval (OSP)
Special Seismic Certification is preapproved under OSP OSP Number:
OR CODE C
Certification Method
Testing in accordance with: ICC-ES AC156 FM 1950-16
Other(s) (Please Specify): OPM-0219
*Use of criteria other than those adopted by the California Building Standards Code, 2022 (CBSC 2022) for component supports and attachments are not permitted. For distribution system, interior partition wall, and suspended ceiling seismic bracings, test criteria other than those adopted in the CBSC 2022 may be used when approved by HCAI prior to testing.
X Analysis
Experience Data DATE: 11/25/2023
Combination of Testing, Analysis, and/or Experience Data (Please Specify):
OPVIA DE CODE
HCAI Approval
Date: 11/25/2023
Name: William Staehlin Title: Senior Structural Engineer
Condition of Approval (if applicable):

"A healthier California where all receive equitable, affordable, and quality health care"

HCAi

STATE OF CALIFORNIA - HEALTH AND HUMAN SERVICES AGENCY



EQUIPMENT ANCHORAGE & SEISMIC ENGINEERING

5877 Pine Ave, Ste. 210 Chino Hills, CA. 91709 Phn: (909) 606-7622

The Department of Health Care Access and Information
PREAPPROVAL OF MANUFACTURER'S CERTIFICATION
OPM-0219

THIS PREAPPROVAL CONFORMS TO THE 2022 CALIFORNIA BUILDING CODE

MANUFACTURER:

BELIMED INFECTION CONTROL

EQUIPMENT NAME: ULTRASONIC WASHERS


Sheet: 1 of 8 Date: 9/18/23

GENERAL NOTES

- 1. THIS HCAI PREAPPROVAL OF MANUFACTURER'S CERTIFICATION (OPM) IS BASED ON THE 2022 CBC. THE DEMANDS (DESIGN FORCES) FOR USE WITH THIS OPM SHALL BE BASED ON THE 2022 CBC
- 2. THIS DOCUMENT MAY ONLY BE USED WITH THE EXPRESS WRITTEN CONSENT OF THE MANUFACTURER LISTED ABOVE FOR THE SPECIFIC PROJECT SITE AND INSTALLATION LOCATION. THIS DOCUMENT IS INVALID WITHOUT SUCH CONSENT.
- 3. THIS PREAPPROVAL CONFORMS TO THE 2022 CALIFORNIA BUILDING CODE WHERE SDS IS NOT GREATER THAN 2.30.
- 4. FORCES PER ASCE 7-16 SECTION 13.3.1, EQUATIONS 13.3-1, 13.3-2 & 13.3-3, WHERE SDS = 2.30, \mathbf{a}_p = 1.0, \mathbf{I}_p = 1.5, \mathbf{r}_p = 1.5,
- 5. THIS PREAPPROVAL COVERS ONLY THE SUPPORTS AND ATTACHMENTS OF THE EQUIPMENT TO THE STRUCTURE.
- 6. ALL DESIGN FORCES SHOWN ON THE DRAWINGS ARE FACTORED LOADS THAT SHALL BE USED FOR STRENGTH DESIGN.
- 7. CONCRETE SLAB ON METAL DECK DETAIL VALID FOR DEMANDS SHOWN AT ANY ELEVATION IN THE BUILDING. (i.e. z/h < 1)
- 8. CONCRETE SLAB DETAIL VALID FOR DEMANDS SHOWN AT OR BELOW GRADE. (i.e. z/h = 0)

9. RESPONSIBILITIES OF THE STRUCTURAL ENGINEER OF RECORD OF THE BUILDING

- A. PROVIDE SUPPORTING STRUCTURE TO SUPPORT WEIGHTS AND FORCES SHOWN IN ADDITION TO ALL OTHER LOADS.
- B. VERIFY THAT THE INSTALLATION IS IN CONFORMANCE WITH THE 2022 CBC AND WITH THE DETAILS, MATERIAL AND GAGE OF THE UNIT WHERE ATTACHMENTS ARE MADE AGREE WITH THE INFORMATION SHOWN ON THE PREAPPROVAL DOCUMENTS.
- C. VERIFY THAT PROJECT SPECIFIC VALUES OF SDS & z/h RESULT IN SEISMIC FORCES (Eh, Ev) THAT DO NOT EXCEED THE VALUES ON THE DETAILS.
- D. VERIFY THAT THE CONCRETE SLAB TO WHICH THE EQUIPMENT IS ANCHORED MEETS THE REQUIREMENTS OF THE APPLICABLE ICC ESR REPORT. AND THIS OPM.
- E. VERIFY THAT THE ANCHORS ARE AN ADEQUATE DISTANCE FROM ANY SLAB EDGES OR OPENINGS (SEE TYPICAL DETAIL ON SHEET 2).
- F. VERIFY THAT ALL NEW OR EXISTING ANCHORS ARE AN ADEQUATE DISTANCE FROM THE UNIT ATTACHMENTS AND CHECK FOR INTERACTION WHERE OTHER ANCHORS ARE WITHIN 18" OR 6hef FROM THIS UNIT'S ANCHORS.

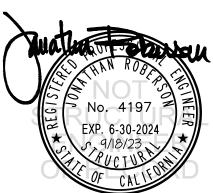

EQUIPMENT ANCHORAGE & SEISMIC ENGINEERING

www.EquipmentAnchorage.com

BELIMED INFECTION CONTROL

ULTRASONIC WASHERS

DES. J. ROBERSON

JOB NO. 11-2319

DATE 9/18/23

2 2

8 sheets


10. EXPANSION ANCHORS:

A. ATTACHMENT IS TO BE MADE WITH THE ANCHORS LISTED BELOW AND INSTALLED AS DESCRIBED IN THE CORRESPONDING ICC REPORT.

Anchor Diamete		Min. f'c (psi)	Anchor Type	ICC Report No.	Min. Embed.	Min. Spacing	Min. Edge Dist.	Min. Conc. Thickness	Torque Test	Direct Tension Test
3/8"	Sand Light Weight	3000	Hilti Kwik Bolt TZ2	ESR-4266	2"	6.75"	12"	See Detail "A"	30 FT-LB	N/A
1/2"	Normal Weight	3000	Hilti Kwik Bolt TZ2	ESR-4266	2"	8"	28	4"	50 FT-LB	1605 lb

- B. THIS PREAPPROVAL ALLOWS FOR UP TO A MAXIMUM OF 2 ADJACENT CONCRETE SLAB EDGES, 28" AWAY MINIMUM (i.e. CORNER).

 SEE ADJACENT DETAIL FOR ADDITIONAL MINIMUM ALLOWABLE CONCRETE EDGE DISTANCES.
- C. TESTING AND SPECIAL INSPECTION OF EXPANSION ANCHORS SHALL BE PERFORMED BY AN APPROVED INDEPENDENT AGENCY EMPLOYED BY THE FACILITY OWNER PER CBC 1704A & 1910A.5 AND CAC 7-149. ALL REPORTS SHALL BE SENT TO THE INSPECTOR OF RECORD, OWNER AND THE ARCHITECT OR ENGINEER IN RESPONSIBLE CHARGE.
 - (i) AFTER AT LEAST 24 HOURS HAVE ELAPSED SINCE INSTALLATION, DIRECT PULL TENSION TEST OR TORQUE TEST AT LEAST 50% OF THE ANCHORS.
 - (ii) ACCEPTANCE CRITERIA:
 - DIRECT TENSION TEST: THE ANCHOR SHOULD HAVE NO OBSERVABLE MOVEMENT AT THE TEST LOAD. A PRACTICAL WAY TO DETERMINE OBSERVABLE MOVEMENT IS THAT THE WASHER BECOMES LOOSE.
 - TORQUE TEST: THE APPLICABLE TORQUE MUST BE ACHIEVED WITHIN THE FOLLOWING LIMITS: WEDGE TYPE: 1/2 TURN OF THE NUT
 - (iii) IF ANY ANCHOR FAILS, TEST ALL ANCHORS.
- D. AVOID DAMAGING EXISTING STEEL REINFORCING IN CONCRETE SLAB WHEN INSTALLING CONCRETE EXPANSION ANCHORS.
- E. PROVIDE FOR FULL THREAD ENGAGEMENT OF NUT & WASHER.
- 11. BOLTS THROUGH CONCRETE ON METAL DECK
 - A. BOLTS SHALL BE TORQUED BY 3/4 TURN OF THE NUTS AFTER THE SNUG TIGHT (THE SNUG-TIGHT CONDITION IS DEFINED AS THE TIGHTNESS REQUIRED TO BRING THE CONNECTED PLIES INTO FIRM CONTACT) CONDITION IS ACHIEVED, UNLESS OTHERWISE NOTED.
 - B. THROUGH BOLT HOLES SHALL BE 1/16" LARGER THAN BOLT SIZE (HOLE SIZE = BOLT SIZE + 1/16) FOR CONCRETE.
 - C. THROUGH-BOLTS IN CONCRETE SHALL RECEIVE SPECIAL INSPECTION AND TESTING (THROUGH BOLTS WITH STEEL TO STEEL CONNECTION IN TENSION DO NOT REQUIRE TENSION TESTING) IN ACCORDANCE WITH REQUIREMENTS FOR POST-INSTALLED ANCHORS.

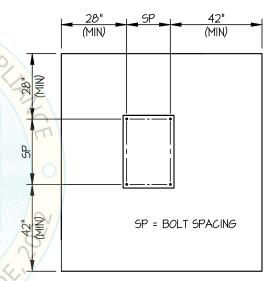

TYPICAL CONCRETE EDGE DETAIL (SLAB ON GRADE ONLY)



EQUIPMENT ANCHORAGE & SEISMIC ENGINEERING

www.EquipmentAnchorage.com

BELIMED INFECTION CONTROL

ULTRASONIC WASHERS MODELS 200c & UW565 DES. J. ROBERSON

JOB NO. 11-2319


DATE 9/18/23

3

8 знеетз

SEISMIC ANCHORAGE

CONCRETE SLAB

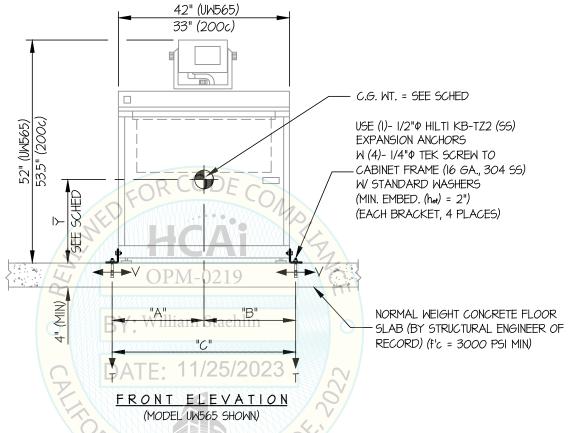

NOTES:


1. FORCES ARE DETERMINED PER 2022 CALIFORNIA BUILDING CODE AND ASCE 7-16. STRENGTH DESIGN IS USED. (EXAMPLE: SDS = 2.30, Δp = 1.0, |p| = 1.5, Rp = 1.5, Ω_0 = 2.0, z/h = 0)

HORIZONTAL FORCE (En) = 1.035 Wp

HORIZONTAL FORCE (Emh) = 2.07 Wp (FOR CONCRETE ANCHORAGE)

VERTICAL FORCE (Ev) = 0.46 Wp

- 2. THIS PREAPPROVAL ENCOMPASSES WEIGHTS AND VERTICAL CG POSITIONS UP TO THE VALUES SHOWN.
- 3. THIS PREAPPROVAL WAS PREPARED WITHOUT KNOWLEDGE OF ANY SITE CONDITION. COMPATIBILITY FOR USE WITH A SITE SHALL BE EVALUATED BY THE STRUCTURAL ENGINEER OF RECORD OF THE INSTALLATION (SEOR). USE REQUIRES APPROVAL BY THE SEOR.
- 4. STRUCTURAL ENGINEER OF RECORD FOR THE INSTALLATION SHALL VERIFY ALL CONDITIONS, EVALUATE INTERACTION WITH ADJACENT EQUIPMENT AND ANCHORS, AND PROVIDE SUPPORT STRUCTURE DESIGNED TO SUPPORT WEIGHTS AND FORCES SHOWN IN COMBINATION WITH ALL OTHER LOADS THAT MAY BE PRESENT.

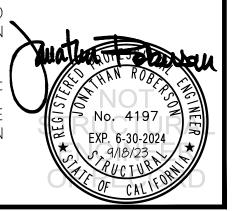

EQUIPMENT ANCHORAGE & SEISMIC ENGINEERING

www.EquipmentAnchorage.com

BELIMED INFECTION CONTROL

ULTRASONIC WASHERS MODELS 200c & UW565 DES. J. ROBERSON

JOB NO. 11-2319

DATE

9/18/23

4


SHEET

8 знеетз

SEISMIC ANCHORAGE

CONCRETE SLAB

USE (I)- I/2" PHILTI KB-TZ2 (SS)
EXPANSION ANCHORS
W (4)- I/4" PTEK SCREW TO
CABINET FRAME (I6 GA., 304 SS):
W/ STANDARD WASHERS
(MIN. EMBED. (Net) = 2")
(EACH BRACKET, 4 PLACES)

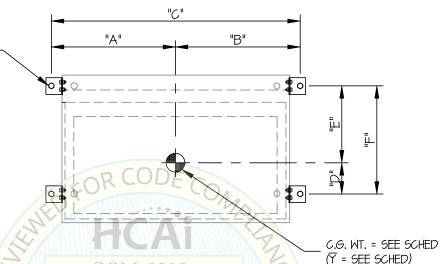

(MODEL UM565 SHOWN)

BY: William Staehlin

MODEL	WEI <mark>GHT (lb.)</mark> MAX	₹ (in.)	"A" (in.)	"B" (in.)	"C" (in.)	"D" (in.)	"E" (in.)	"F" (in.)	++ Tu (lb.)	Yu (lb.)
200c	550	20.7	18.5	18.5	37	6	/ 14	20	638	484
* UW565	855	21.4	23	23	46	5.8	14.2	20	989	761

^{*} THIS UNIT IS USED IN CALCULATION



⁺⁺ VALUES INCLUDE Ω₀

EQUIPMENT ANCHORAGE & SEISMIC ENGINEERING

www.EquipmentAnchorage.com

BELIMED INFECTION CONTROL

ULTRASONIC WASHERS MODELS 200c & UW565 DES. J. ROBERSON

JOB NO. 11-2319

DATE

9/18/23


5

SHEET

of 8 sheets

SEISMIC ANCHORAGE

CONCRETE SLAB ON METAL DECK

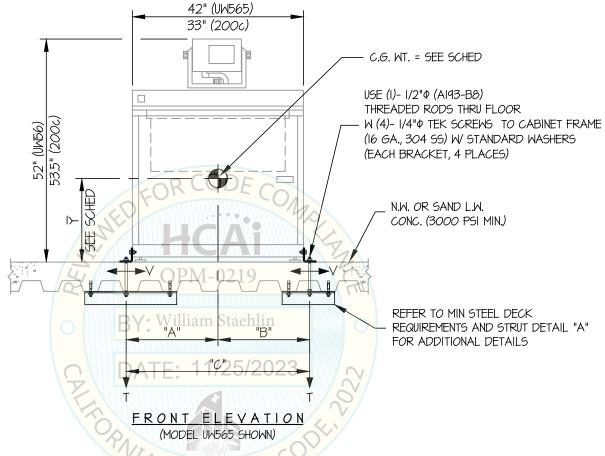

NOTES:


1. FORCES ARE DETERMINED PER 2022 CALIFORNIA BUILDING CODE AND ASCE 7-16. STRENGTH DESIGN IS USED. (EXAMPLE: SDS = 2.30, 20 = 1.0, 20 = 1.5, 20 = 2.0, 20 = 2.0, 20 = 1.5, 20 = 2.0, 20

HORIZONTAL FORCE (En) = 2.76 Wp

HORIZONTAL FORCE (Emh) = 5.52 Wp (FOR CONCRETE ANCHORAGE)

VERTICAL FORCE (Ev) = 0.46 Wp

- 2. THIS PREAPPROVAL ENCOMPASSES WEIGHTS AND VERTICAL CG POSITIONS UP TO THE VALUES SHOWN.
- 3. THIS PREAPPROVAL WAS PREPARED WITHOUT KNOWLEDGE OF ANY SITE CONDITION. COMPATIBILITY FOR USE WITH A SITE SHALL BE EVALUATED BY THE STRUCTURAL ENGINEER OF RECORD OF THE INSTALLATION (SEOR), USE REQUIRES APPROVAL BY THE SEOR.
- 4. STRUCTURAL ENGINEER OF RECORD FOR THE INSTALLATION SHALL VERIFY ALL CONDITIONS, EVALUATE INTERACTION WITH ADJACENT EQUIPMENT AND ANCHORS, AND PROVIDE SUPPORT STRUCTURE DESIGNED TO SUPPORT WEIGHTS AND FORCES SHOWN IN COMBINATION WITH ALL OTHER LOADS THAT MAY BE PRESENT.

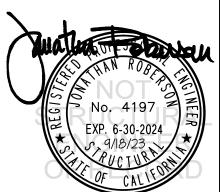

EQUIPMENT ANCHORAGE & SEISMIC ENGINEERING

www.EquipmentAnchorage.com

BELIMED INFECTION CONTROL

ULTRASONIC WASHERS <u> MODELS 200c & UW565</u> DES. J. ROBERSON

11-2319 JOB NO.

9/18/23 DATE


SHEET

SHEETS

SEISMIC ANCHORAGE

CONCRETE SLAB ON METAL DECK

USE (I)- 1/2"Φ (A193-B8) THREADED RODS THRU FLOOR W (4)- I/4" TEK SCREWS TO CABINET FRAME -(16 GA., 304 SS) W/ STANDARD WASHERS (EACH BRACKET, 4 PLACES)

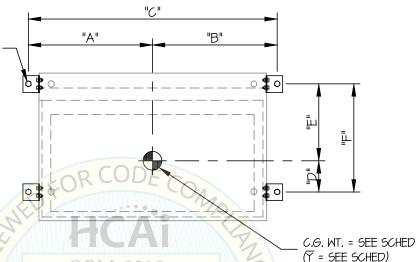

PLAN 2AT BASE (MODEL UM565 SHOWN)

BY: William Staehlin

MODEL	WE <mark>IGHT (</mark> lb.) (MAX)	▼ (in.)	"A" (in.)	"B" (in.)	"C" (in.)	"D" (in.)	"E" (in.)	"F" (in.)	++ Tu (lb.)	++ Vu (lb.)
200c	550	20.7	18.5	18.5	37	6	14	20	879	645
* UW565	855	21.4	23	23	46	5.8	14.2	20	1363	1015

THIS UNIT IS USED IN CALCULATION



11/25/2023

⁺⁺ VALUES DO NOT INCLUDE $\Omega_{
m o}$

EQUIPMENT ANCHORAGE & SEISMIC ENGINEERING

www.EquipmentAnchorage.com

OF

BELIMED INFECTION CONTROL

ULTRASONIC WASHERS MODELS 200c & UW565 DES. J. ROBERSON

JOB NO. 11-2319


DATE 9/18/23

7

8 SHEETS

SEISMIC SUPPORTS & ATTACHMENTS

CONCRETE DETAIL

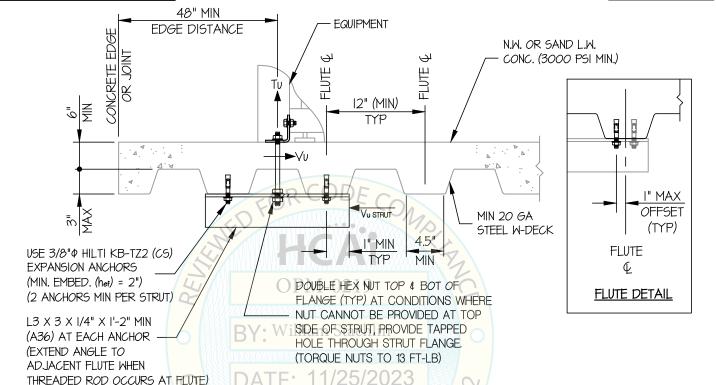

MIN STEEL DECK REQUIREMENTS AND STRUT DETAIL.

PNIA BUIL



EQUIPMENT ANCHORAGE & SEISMIC ENGINEERING

www.EquipmentAnchorage.com

BELIMED INFECTION CONTROL

ULTRASONIC WASHERS MODELS 200c & UW565

DES. J. ROBERSON


JOB NO. 11-2319

DATE 9/18/23

8HEET

ог 8 знеетз

BRACKET DETAIL

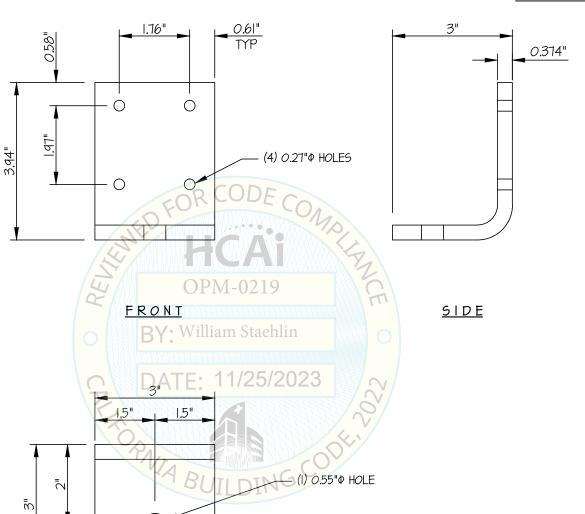

304 S.S. (FY= 35 KSI)

<u>PLAN</u>

